


1. Propyl gallate
2. THBP
3. TBHQ
4. NDGA
5. BHA
6. Ionox 100
7. Octyl gallate
8. BHT
9. Dodecyl gallate

Gradient Program		
Time (min)	%A	%B
0	70	30
10	0	100
20	70	30

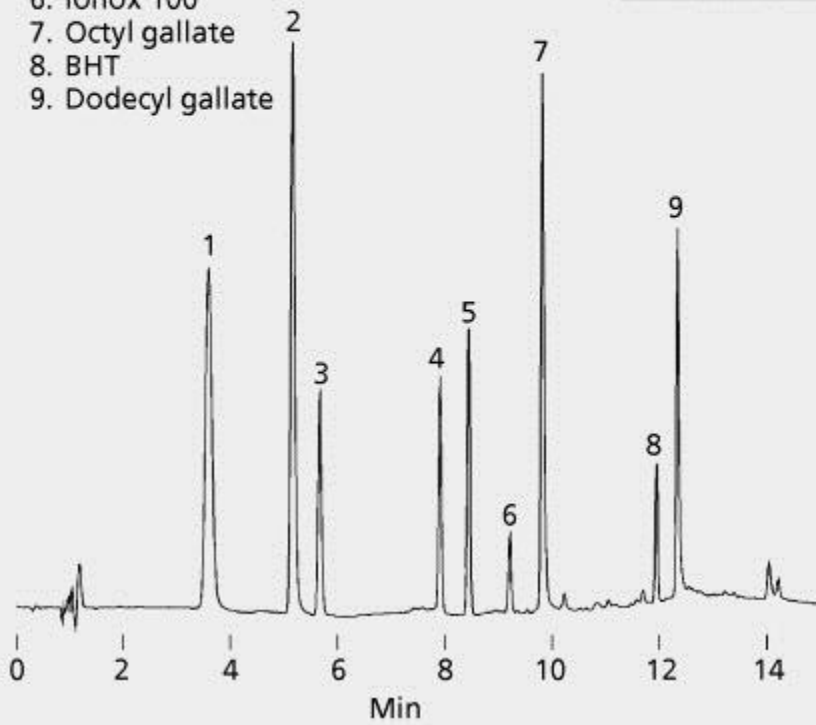

test parameters	column: SUPELCOSIL LC-18, 15 cm × 4.6 mm I.D., 5 µm particles (58230-U)
	mobile phase: (A) 5% acetic acid in deionized water; (B) acetonitrile:methanol (1:1) 70% A/30% B to 100% B, linear gradient over 10 min, hold 10min
	flow rate: 2 mL/min
	detector: UV, 280 nm
	injection: 10 µL, 20 µg/mL each antioxidant